

How to display the Job field (Assay ID) in a TargetLynx dataset

1. Choose File >Report Format

The screenshot shows the TargetLynx XS interface. The 'File' menu is open, and 'Report Format...' is highlighted. The background shows a data table for 'NIC2 (1)' with columns: Std. Conc, RT, Area, IS Area, Response, Primar..., pg/uL, %Dev, and Conditions. Below the table is a chromatogram plot with a red peak labeled 'MRM of 8 channels, ES+ 163.134 > 131.97 1.771e+007'.

Std. Conc	RT	Area	IS Area	Response	Primar...	pg/uL	%Dev	Conditions
1.000	5.10	28910.857		28910.857	bdl			
2.000	5.10	2151542.750		2151542.750	bb	4.3	115.7	
4.000	5.08	2039290.250		2039290.250	bb	4.1	2.1	
8.000	5.08	3278734.500		3278734.500	bb	6.6	-17.4	
	5.01	222.756		222.756	bbl			
	4.98	154.536		154.536	bbl			
	5.04	129.101		129.101	bbl			
	5.08	50.804		50.804	bbl			
	5.07	126.833		126.833	bbl			
	5.05	101.009		101.009	bbl			
	5.12	77.548		77.548	bbl			
	4.85	136.598		136.598	bbl			
	5.00	59.717		59.717	bbl			

2. Enable Experiment Report and select 'Print Selection for All Samples' and 'Header':

The 'Report Format Options' dialog box is shown. The 'Experiment Report' checkbox is checked. Under 'Print Only', 'Print Selection For All Samples' is selected. Under 'Options', 'Header' is checked.

Report Format Options

Compounds Report	Samples Report	Method Report	Totals Report	Audit Report
Statistics Report	Sample Template Report	Overview Report	PICS Report	Matrix Factor Report
General	Compound Summary Report	Sample Summary Report	Calibration Report	Experiment Report

Enable

Orientation

Portrait
 Landscape

Header Fields...

Print Only


Print Full Record For All Samples
 Print Full Record For First Sample Only
 Print Selection For All Samples
 Print Selection For First Sample Only
 Print Full Record For First Sample And Selection For Rest

Options

Pump
 Header
 Autosampler
 Calibration
 Detector(s)
 MS
 GC
 Function


OK Cancel Help

3. Click Header Fields and add the Job Code field:

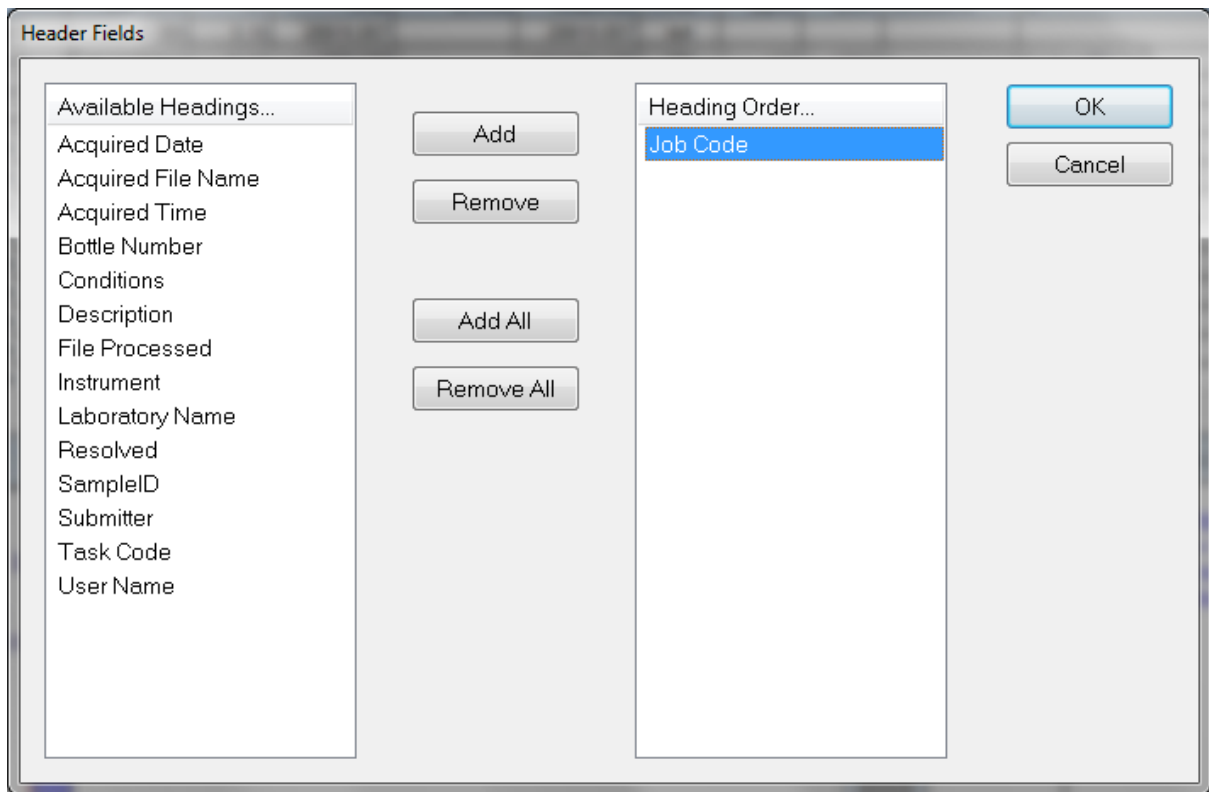

4. Click OK (twice) to close the dialogs.

5. Choose File > Print Preview

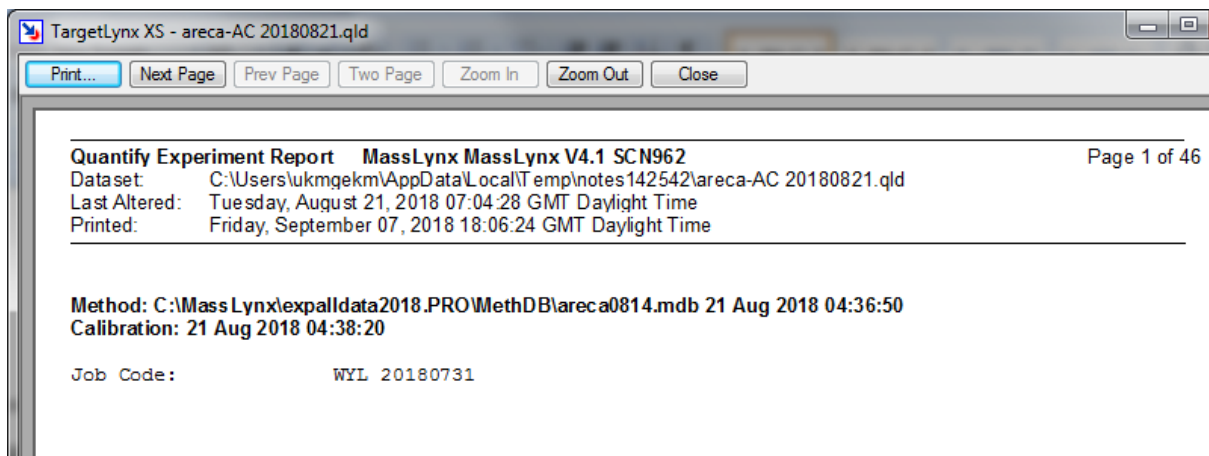

6. Click 'Next Page' to scroll through samples. To use the LIMS Interface to send the results to LIMS or Middleware, the Job Code must be the same for all samples.

Note: if there are two or more different Job Codes, it will not be possible to upload the results.